



## HiCrome™ Bacillus Agar

M1651

### Intended use

Recommended for isolation and differentiation between various species of *Bacillus* from a mixed culture in foods, clinical and non-clinical samples by chromogenic method.

### Composition\*\*

Ingredients	Gms / Litre
Peptone	10.000
HM extract #	1.000
D-Mannitol	10.000
Sodium chloride	10.000
Chromogenic mixture	3.200
Phenol red	0.025
Agar	15.000
Final pH ( at 25°C)	7.1±0.2

\*\*Formula adjusted, standardized to suit performance parameters

#-Equivalent to Meat extract

### Directions

Suspend 49.22 grams in 1000 ml purified/distilled water. Heat to boiling to dissolve the medium completely. Sterilize by autoclaving at 15 lbs pressure (121°C) for 15 minutes. Cool to 50°C and aseptically add rehydrated contents of 1 vial of Bacillus Selective Supplement (FD324) if desired. Mix well and pour into sterile Petri plates.

### Principle And Interpretation

Majority of *Bacillus* species apparently have little or no pathogenic potential and are rarely associated with disease in humans or lower animals. The principal exception to this are *Bacillus anthracis*, the agent of anthrax, and *Bacillus cereus*, but a number of other species, particularly those of the *B.subtilis* group, have been implicated in food poisoning and other human and animal infections (7). *Bacillus cereus* causes food poisoning due to consumption of contaminated rice (2,5,10), other starchy foods such as potato, pasta and cheese have also been implicated, eye infections and a wide range of other clinical conditions like abscess formation, meningitis, septicemia and wound infection.

HiCrome™ Bacillus Agar is based on the formulation of MYP Agar formulated by Mossel et al (6) used for enumeration of *Bacillus cereus* and *Bacillus thuringiensis* when present in large number in certain foodstuffs

The medium contains peptone and HM extract, which provide nitrogenous, carbonaceous compounds, long chain amino acids, vitamins and other essential growth nutrients. Mannitol serves as the fermentable carbohydrate, fermentation of which can be detected by phenol red. Mannitol fermenting organisms like *B. megaterium* yield yellow coloured colonies. The chromogenic mixture present in the medium is cleaved by the enzyme beta-glucosidase found in *B.cereus* resulting in the formation of blue colonies. *B.thuringiensis* also grows as blue/green colonies on this medium as *B.cereus* and *B.thuringiensis* are biochemically identical, however *B.cereus* shows flat colonies with distinct blue centers, while *B.thuringiensis* shows irregular margins. If selective isolation of *B.cereus* or *B.thuringiensis* is required aseptically add Bacillus Selective supplement (FD324).

### Type of specimen

Clinical samples - Blood, pus, eye exudates; Food and dairy samples.

### Specimen Collection and Handling:

For clinical samples follow appropriate techniques for handling specimens as per established guidelines (3,4).

For food and dairy samples, follow appropriate techniques for sample collection and processing as per guidelines (1,8,9).

After use, contaminated materials must be sterilized by autoclaving before discarding.

## Warning and Precautions :

In Vitro diagnostic Use only. Read the label before opening the container. Wear protective gloves/protective clothing/eye protection/ face protection. Follow good microbiological lab practices while handling specimens and culture. Standard precautions as per established guidelines should be followed while handling clinical specimens. Safety guidelines may be referred in individual safety data sheets.

## Limitations :

1. Due to variable nutritional requirements, some strains may show poor growth on this medium.
2. Slight colour variation may be observed depending upon the utilization of the substrate by the organism.

## Performance and Evaluation

Performance of the medium is expected when used as per the direction on the label within the expiry period when stored at recommended temperature.

## Quality Control

### Appearance

Cream to yellow homogeneous free flowing powder

### Gelling

Firm, comparable with 1.5% Agar gel

### Colour and Clarity of prepared medium

Red coloured, clear to slightly opalescent gel forms in Petri plates

### Reaction

Reaction of 4.92% w/v aqueous solution at 25°C. pH : 7.1±0.2

### pH

6.90-7.30

### Cultural Response

Cultural characteristics observed after an incubation at 30°C for 24-48 hours.

### Cultural Response

Organism	Inoculum (CFU)	Growth w/o addition of FD324	Recovery w/ o addition of FD324	Growth w/ addition of FD324	Recovery w/ addition of FD324	Colour of Colony
<b>Cultural Response</b>						
<i>Bacillus subtilis subsp. spizizenii</i> ATCC 6633 (00003*)	50-100	fair	20-30%	inhibited	0%	yellowish green to green colonies
<i>Bacillus cereus</i> ATCC 10876	50-100	good-luxuriant	>=50%	good-luxuriant	>=50%	light blue, large, flat colonies with blue centre
<i>Bacillus thuringiensis</i> ATCC 10792	50-100	good-luxuriant	>=50%	good-luxuriant	>=50%	light blue, large, flat colonies with irregular margins
<i>Bacillus megaterium</i> ATCC 14581	50-100	good-luxuriant	>=50%	inhibited	0%	yellow, mucoid colonies
<i>Bacillus coagulans</i> ATCC 7050 (00002*)	50-100	good-luxuriant	>=50%	inhibited	0%	pink, small, raised colonies
<i>Bacillus pumilus</i> ATCC 14884	50-100	good-luxuriant	>=50%	poor	10-20%	light green to green colonies
<i>Staphylococcus aureus subsp. aureus</i> ATCC 25923 (00034*)	50-100	luxuriant	>=50%	inhibited	0%	yellow colonies
<i>Enterococcus faecalis</i> ATCC 29212 (00087*)	50-100	luxuriant	>=50%	inhibited	0%	light green to green colonies

\* - Corresponding WDCM numbers

Please refer disclaimer Overleaf.

## Storage and Shelf Life

Store dehydrated medium in a tightly closed container and the prepared medium at 2-8°C. Use before expiry date on the label. On opening, product should be properly stored dry, after tightly capping the bottle in order to prevent lump formation due to the hygroscopic nature of the product. Improper storage of the product may lead to lump formation. Store in dry ventilated area protected from extremes of temperature and sources of ignition. Seal the container tightly after use. Use before expiry date on the label.

Product performance is best if used within stated expiry period.

## Disposal

User must ensure safe disposal by autoclaving and/or incineration of used or unusable preparations of this product. Follow established laboratory procedures in disposing of infectious materials and material that comes into contact with clinical sample must be decontaminated and disposed of in accordance with current laboratory techniques (3,4).

## Reference

1. American Public Health Association, Standard Methods for the Examination of Dairy Products, 1978, 14th Ed., Washington D.C.
2. Bouza E., Grant S., Jordan C. et al, 1979, Arch. Ophthalmol., 97:488.
3. Isenberg, H.D. Clinical Microbiology Procedures Handbook. 2<sup>nd</sup> Edition.
4. Jorgensen, J.H., Pfaller, M.A., Carroll, K.C., Funke, G., Landry, M.L., Richter, S.S and Warnock, D.W. (2015) Manual of Clinical Microbiology, 11th Edition. Vol. 1.
5. Mortimer P. R. and McCann G., 1974, Lancet, 1043.
6. Mossel D. A. A., Koopman M. J. and Jongerium E., 1967, Appl. Microbiol., 15:650.
7. Murray P. R., Baron J. H., Pfaller M. A., Jorgensen J. H. and Tenover F. C., (Eds.), 2003, Manual of Clinical Microbiology, 8th Ed. American Society for Microbiology, Washington, D.C.
8. Salfinger Y., and Tortorello M.L. Fifth (Ed.), 2001, Compendium of Methods for the Microbiological Examination of Foods, 5th Ed., American Public Health Association, Washington, D.C.
9. Wehr H. M. and Frank J. H., 2004, Standard Methods for the Microbiological Examination of Dairy Products, 17th Ed., APHA Inc., Washington, D.C.
10. Wohlgenuth K., Kirkbride C. A., Bicknell E. J. and Ellis R. P., 1972 Am. Vet. Met. Ass. 161:1691.

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
Revision : 03 / 2017



In vitro diagnostic medical device



CE Marking



Storage temperature



Do not use if package is damaged



HiMedia Laboratories Pvt. Limited,  
23 Vadhani Industrial Estate,  
LBS Marg, Mumbai-86, MS, India



CE Partner 4U, Esdoornlaan 13, 3951  
DB Maarn The Netherlands,  
[www.cepartner4u.eu](http://www.cepartner4u.eu)

Please refer disclaimer Overleaf.

**Disclaimer :**

User must ensure suitability of the product(s) in their application prior to use. Products conform solely to the information contained in this and other related HiMedia™ publications. The information contained in this publication is based on our research and development work and is to the best of our knowledge true and accurate. HiMedia™ Laboratories Pvt Ltd reserves the right to make changes to specifications and information related to the products at any time. Products are not intended for human or animal or therapeutic use but for laboratory, diagnostic, research or further manufacturing use only, unless otherwise specified. Statements contained herein should not be considered as a warranty of any kind, expressed or implied, and no liability is accepted for infringement of any patents.